



Quick Start Guide


StrataVac
FOR 1 HEAD SHORT PATH KITS

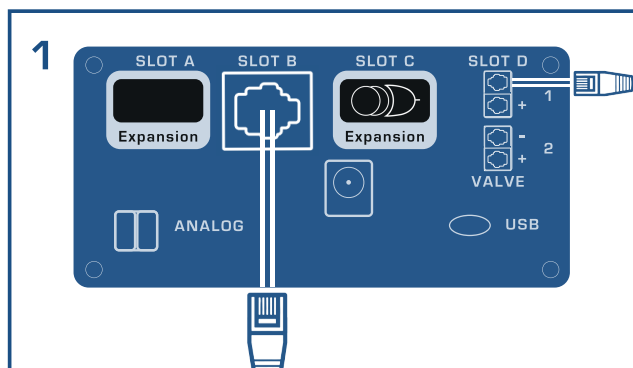


1 Connect sensor cable to Slot B and connect the valve to Slot D (top)

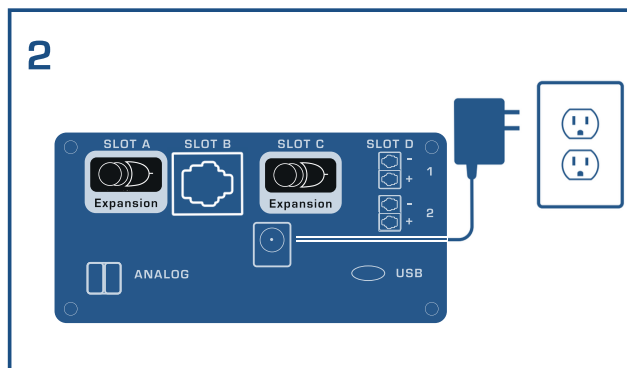


2 Plug StrataVac into wall outlet using the power supply supplied by DigiVac

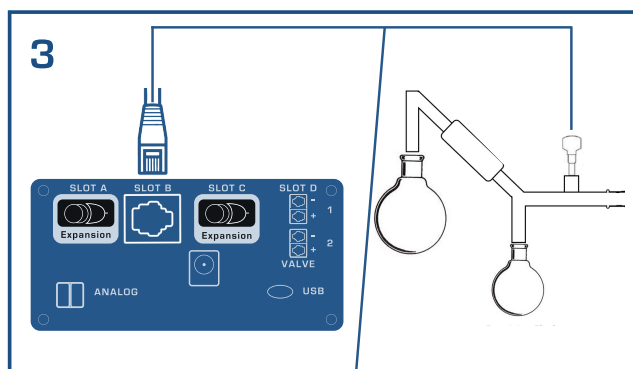


3 Take sensor and mount it as close to the reaction as possible

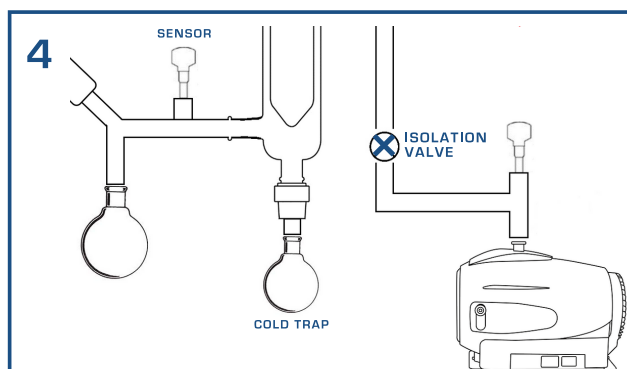

4 Place your isolation valve between the cold trap and your pump

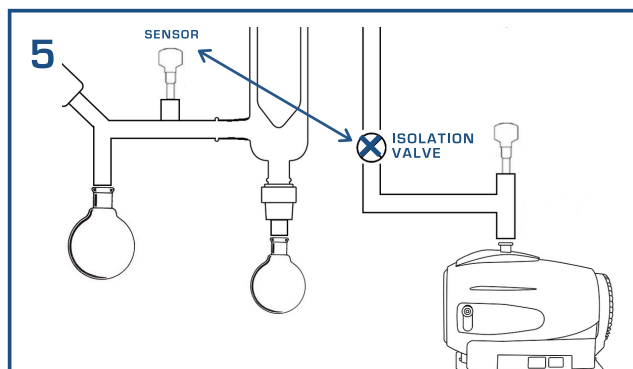

5 The valve will open and close based on the pressures that Sensor B reads

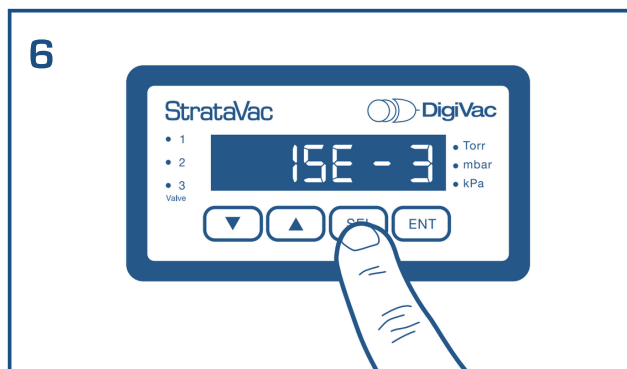

6 You are ready to measure and control with the StrataVac!

Quick Start Guide

StrataVac



To navigate to the upper set point, press Select (SEL) until you see “U” on the display.



Use the up/down arrows to change the set point. Once the desired set point is reached, press Enter (ENT) to save.



To navigate to the lower set point, press Select (SEL) until you see “L” on the display.



Repeat step two to set and save the lower set point.